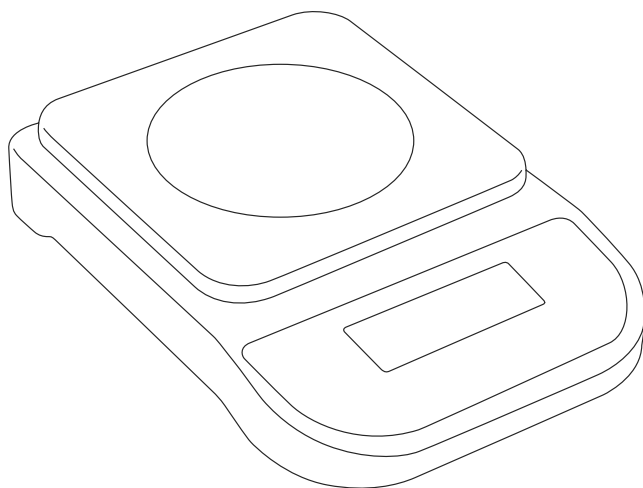
GRAM

SERIES / SERIE / SÉRIE

SPX

600D / 1000D / 3000D / 6000D

EN

ES

FR

CE

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Installation	5
Keyboard description	6
Symbols on display	7
Operations	8
Turn on the scale	8
Reset to zero	8
Tare function	8
Weight unit selection	8
Pieces counting function	9
Turn off and automatic turn off	9
Calibration	10
Guarantee	11

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ENGLISH

INSTALLATION

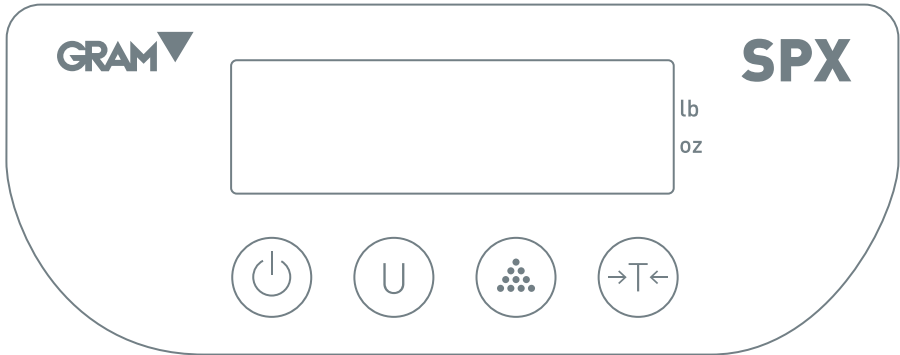
Recommendations for maximum optimization.

To obtain both the best product possible and the best results, we strongly recommend installing the scale by following these guidelines below:

- Install the scale on a flat surface.
- Place the scale on a stable and vibration free surface.
- Do not expose to direct sunlight
- Do not use the scale in an atmosphere containing combustible corrosive gas.
- Do not store it in a dusty or dirty place.
- Store at room temperature 59° - 86° F (15° - 30° C)
- Use in a room with less than 95% RH humidity
- Keep the scale away from objects that may cause magnetic or acoustic interference.
- Calibrate it in its final location and verify from time to time it did not move
- because of the optional calibration weighing.

VERY IMPORTANT: Keep the original package. It could be useful in the future in transporting the balance.

KEYBOARD DESCRIPTION

To turn on or turn off the balance.

To select the weight unit.

Press it to enter to counting mode.

Reset the scale to zero or tare.

SYMBOLS ON DISPLAY

— Negative value.

▶ Units indicator.

g Grams units.

pcs Counting mode.

Zero Zero indicator.

○ Stable.

OPERATIONS

Turn on the scale

Press the  key to turn on the scale. The scale will show « 88888 », and when it will be ready for use.

Reset to zero

If the weighing pan is empty and the reading is not « 0 » or the zero indicator isn't displayed, press the  key to display « 0 » and the zero indicator appears.

Tare function

This function allows obtaining the net weight of a product.

1. With display showing zero value, place an empty container on the weighing pan.
2. Press the  key.
3. The balance will subtract the container's weight, put the product into the container, after stable, the value displayed is the weight of the product (net weight).

To return to normal weighing mode, remove the container from platter and press the  key.

Weight unit selection

Press the  key to select the desired unit weight.

Pieces counting function

1. Press the  key to turn on the balance.
2. Place on the weighing pan a quantity of items as initial sample.
3. Press and hold the  key until to view “**S-10**”.
4. Press again the  key to select the desired sample number (**S-10, S-20 or S-50**).
5. The balance begin to count after a few seconds.
6. Press the  key to exit.

The greater quantity of pieces put in the initial sample, the greater accuracy in balance's results.

Turn-off and automatic turn-off

To turn-off the balance, press the  key.

If your scale have the auto-power off function, it will automatically turn off after 2 minutes without operation.

Calibration

1. Press the  to turn on the scale for warm-up for 1-2 minutes.
2. Press the  key , the display will show **“0.0”**
3. Press and hold the  key for 3 seconds, the display will show the message **“CAL”** and then the calibration weight value.
4. Place on the weighing pan the calibration weight.
5. The display will show the weight and twinkle and then the message **“PASS”**.
6. The calibration has been completed.

NOTE: If your balance still does not read accurately, it means that the calibration failed and please follow the above steps to re-calibrate it again.

Please, do it on a very stable and flat surface.

GUARANTEE

The scale is warranted against defects in material and workmanship for a period of 1 year from the date of delivery.

During the warranty period, **GRAM PECISION SL** will repair any product that proves to be defective.

Limitation of warranty

This warranty does not apply to defects resulting from misuse of any product or part. This warranty also does not apply to damage from overload or problems arising from failure to follow instructions (Especially concerning the MAINTENANCE TIPS).

The warranty does not cover shipping costs necessary to repair the scale. Therefore, shipping charges to and from the repair location must be paid by the owner.

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