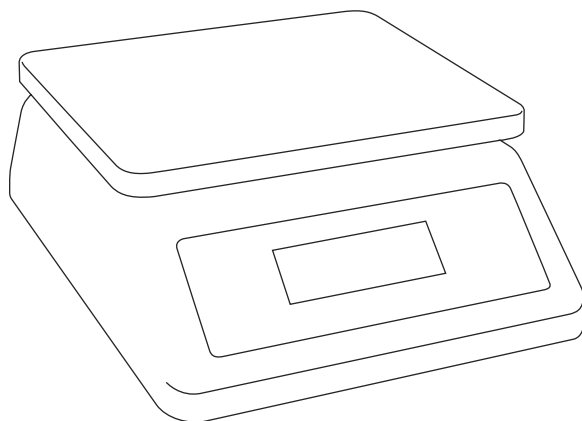
GRAM

SERIES / SERIE / SÉRIE

GX

6000 / 15K / 30K

EN

ES

FR

CE

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Technical characteristics	5
Front keypad	6
backside display	6
Messages	7
Installation	7
Turn-on / Turn-off the balance	8
Reset to zero	8
Tare function	8
Stand-by function	9
State of battery indicators	9
Battery recharging	9
Parameters setup	9
Auto-zero range setting	9
Filter setting	10
Select weighing unit	10
Limits setup	10
Auto-turn-off	11
Buzzer	11
Resolution setup (d)	11
On range setup (on-ra)	12
Displays setup (d-dp)	12
Counting pieces mode	12
Limits indicators	13
Calibration	13
Maintenance	14
Guarantee	15

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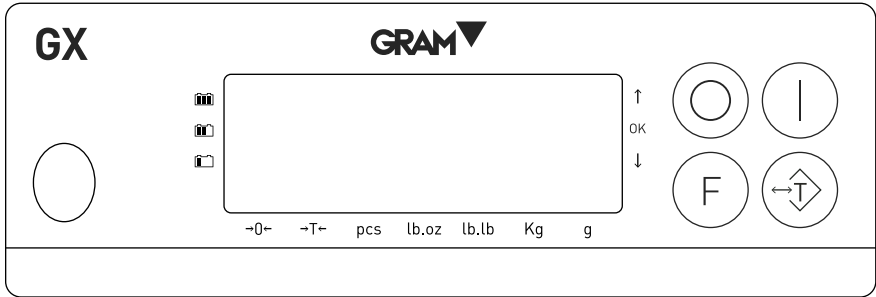
ENGLISH**TECHNICAL CHARACTERISTICS**

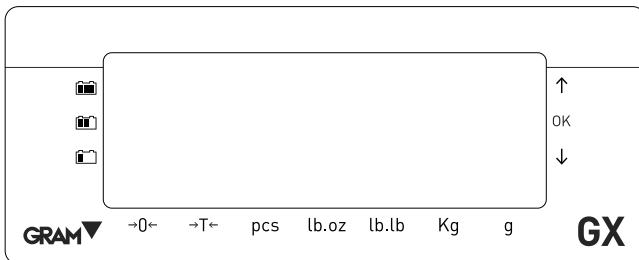
Model	GX-6000	GX-15K	GX-30K
Maximal capacity	6kg	15kg	30kg
Minimal capacity	20e		
Readability	0.5g	1g	2g
Tare range	100% maximum capacity		
Precision	III		
Operating temperature	-5~+35°C		
Storage temperature	-25~+50°C		
Power	6V / 4Ah rechargeable battery and AC/Adaptor		
Dimensions	232 x 284 x 95 mm		
Pan size	225 x 186 mm		
Net weight	3,75 kg		

FRONT KEYPAD

-  To turn on the balance and reset to zero the reading 's display.
-  Tare function, to subtract the weight of a container.
Access to calibration function.
-  To turn off the balance.
-  To setup parameters and data introduction.

BACKSIDE DISPLAY

MESSAGES

- dc x.xx** Shows the current battery voltage “ x.xx V”
- - - - -** It means that the weight placed on the weighing pan exceeds the maximum capacity, the balance will beep, you must immediately remove the weight from the weighing pan.
- AdC-** Means that the digital analogic converter (A/D) values, have been exceeded, the balance will beep.
- Lo-** Indicates that battery voltage is lower than 5.8V. You can go on using the balance for a while , but you should charge it immediately.
- C_End** Indicates that the battery charging is completed.
- Err-1** Battery voltage is lower than 5.6V. The balance will automatically turn offafter this message is shown. Connect the AC/Adaptor and charge it.
- Err-2** Weight adjustment error.

INSTALLATION

Level the balance, turning on the adjustable feet located at the bottom of the scale, so that the bubble remains in the centre of the circle.

TURN ON / TURN OFF THE BALANCE

Turn on

Press the key , the balance will beep. The display will show the version number, the voltage level of battery and it will perform a self-checking 9, 8, 7, ,1,0, finally the display will show the value zero "0", now the balance is ready for weighing.

Turn off.

Keep pressed the key  for few seconds to turn off the balance.

Auto Turn off

After 10 minutes without using the balance, it will automatically turn off.

Turn off for low voltage

When the voltage is less than 5.6V, the display will show the message Err-1 and the balance will turn off.

RESET TO ZERO

When there are no objects placed on the pan and display doesn't indicate "0", press the key , to correct the deviation, now the display reads "0".

TARE FUNCTION

Place a container on the weighing pan, press the key , the display will show zero again and Tare symbol will light on. Put the product inside the

container, the display will show its net weight. When the container is removed, the display will show a negative value and Tare symbol will disappear.

STAND-BY FUNCTION

After 40 seconds of no use, the balance will go into power saving mode (stand-by) and the display will show "0".

STATE OF BATTERY INDICATORS

"High" means that voltage is higher than 6.3V.

"Mid" means that voltage is between 6V and 6.3V.

"Low" means that voltage is lower to 6V.

BATTERY RECHARGING

The internal rechargeable battery is off. When the display shows "-Lo-", turn off the balance and recharge it immediately. The time needed to charge the battery at 100% is 12 hours.

PARAMETERS SETUP

Press the key  for 3 seconds, the display will show "range", continuously pressing the key , will be displayed the following functions: **range \ Units \ A-off \ FILt \ Zero \ Buzzer \ d \ on-ra \ d-dp.**

Auto-zero range setting

Press the  key for 3 seconds until the displays shows the message **"range"**, press again the key  until "2Ero" message is displayed. Press the key , the display will show the current zero range, press the key 

to choose the desired range: 0.5d,1d,1.5d,2d,2.5d,3d,3.5d,4d,4.5d,5d. ("d" is the division value, when n=3000). Press the key  to confirm.

Filter setting

Press the key  for 3 seconds until the message **"range"** is displayed, press again the key  until seeing the message "FILtX" (FILt1 for a stable environment, FILt2 for an unstable environment and FILt3 for very unstable environment).

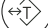
Press the key  to access to three available modes, press the keys  to select the desired mode and confirm pressing key .

Select weighing unit

Press the key  for 3 seconds until the message **"range"** is displayed, press again the key  until the display shows "Unit5", press the key  to access to the parameter, use the key  to choose the desired weighing unit and confirm by pressing .

Available weighing unit: kg, g, Lb.lb, Lb.oz, pcs. (counting-pieces)

Limits setup

Press the key  for 3 seconds until the display shows the message **"range"**, press the key  to enter. The display will show **"off"**, press the key  to select the mode **"on"** and then press . Now the indicator "under" will light on. Press the key  to change the first digit and use to  move to the next digit. Press  to confirm the introduced value.

Now, the indicator "over" will light on. Press the key  to change the first digit and use  to move to the next digit. Press  to confirm the introduced value.

Auto Turn-off (AOFF)

Press the key  for 3 seconds until the display shows the message “range”, press again the key  until the display shows “Aoff”.

Press the key  to enter and press  to select the desired option. To confirm and exit press the key .

N= NO (Auto turn-off disable)

Y= YES (Auto turn-off able)

Buzzer

Press the key  for 3 seconds until the display shows the message “range”, press again  the until you see “Buzzer”

N= NO (Buzzer disabled in over/under menu)

Y=YES (Buzzer activated in over/under menu)

Press the  to enter and press the key  to select the desired option. To confirm and exit press  the key.

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Press the key  for 3 seconds until the display shows the message “range”, press again the key  until you see “d”

Press the  to access to the function and press  to select the desired option. To confirm press the  key.

On range setup (on-ra)

Press the key  for 3 seconds until the display shows the message “range”, press again the key  until you see “on-ra”. Press the  to access to the function  and to select the desired option between 20% of F.E and 100% of F.E. To confirm press the key .

Displays setup (d-dp)

Press the key  for 3 seconds until the screen shows the message “range”, press again the key until you see “d-dp”. Press the key  to access and press the key  to select the desired option.

N= Only one display**Y= Two displays**

Press the key  to confirm and exit.

Counting pieces mode

Enter to “change units” and select the option “pcs”. Press the key  and then the indicator “pcs” will light on.

Sampling procedure:

Press the key , the message “count” will appear. Put the sample on the weighing pan and press the key  in order to introduce the sample’s value. Press the key  to increase the value of the digit and press the key  to move to the right. When the last digit flashes, press the key  to finish the sampling, now the display will show the number of pieces used as a sample. To exit, enter again to “change units” and select the desired unit weight, the

balance will return to normal weighing mode.

LIMITS INDICATORS

- When the weight exceeds the high limit, the indicator “**over**” will light on.
- When the weight is between the low and high limits, the indicator “**accept**” will light on.
- When the weight is below the low limit, the indicator “**low**” will light on.

Note: If the display is instable, the limits indicators won't light on.

Attention: When the low limit is greater than the high limit, the display will show the message Err2, in this case, please cancel the operation.

CALIBRATION

Resolution selection

1. There is a switch on the bottom of the scale protected by a cover and two screws. Move the switch towards the keyboard.
2. Turn on the scale pressing the key .
3. Keep pressed the key  until the display shows the message “rAngE”
4. Press the key  until you see “d”
5. Press the key  to access.
6. Press the key  repeatedly to see the resolution values available.
7. Press the key  to select and confirm.

Adjustment

1. Move the switch in the position closest to keyboard.
2. Press the key  for 3 seconds, the display will show "CAL 0.000"
3. Press the key , the zero adjustment will be completed.
4. The display will show the weight calibration.
5. Place the weight calibration on the weighing pan.
6. Press the  to perform the calibration.
7. Once calibration is completed, move the switch towards the rear part of the balance.

MAINTENANCE

1. If the balance is not to be used for a long period of time, please recharge the battery at least every 3 months.
2. Do not exceed the maximum capacity of the balance.
3. Do not drop objects violently on the weighing pan.
4. The damages caused by overloading are not covered by the warranty.
5. Do not use sharp objects on the keyboard, if the rubber is damaged, the waterproof protection would be damaged.
6. Keep the balance always clean.

GUARANTEE

This balance is guaranteed for one year from the delivery date. The guarantee covers any fabrication defect of the material.

During this period, GRAM PRECISION covers the manpower and the spare parts necessary to repair the scale.

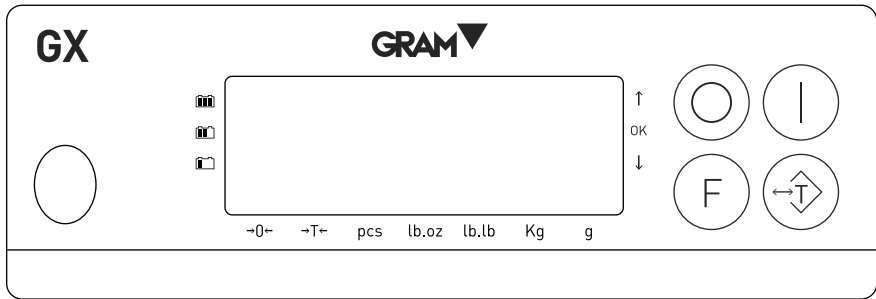
This guarantee does not cover the failures caused by an inappropriate use or overload.

The guarantee does not cover the freight cost (transport) necessary to repair the balance.

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